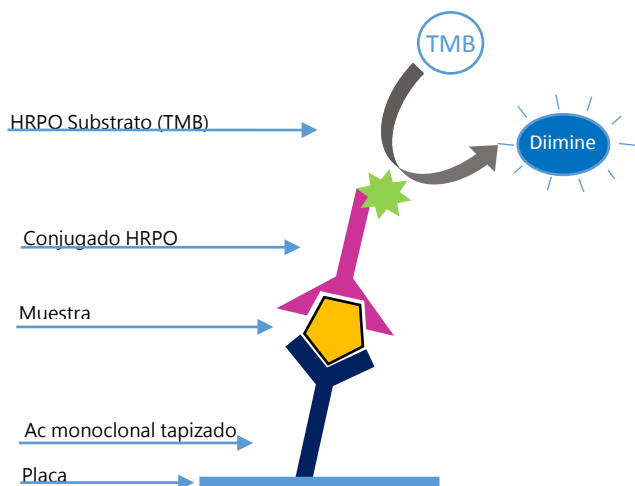
INGEZIM PCV DAS

R.11.CIR.K2

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
INGEZIM PCV DAS

R.11.CIR.K2

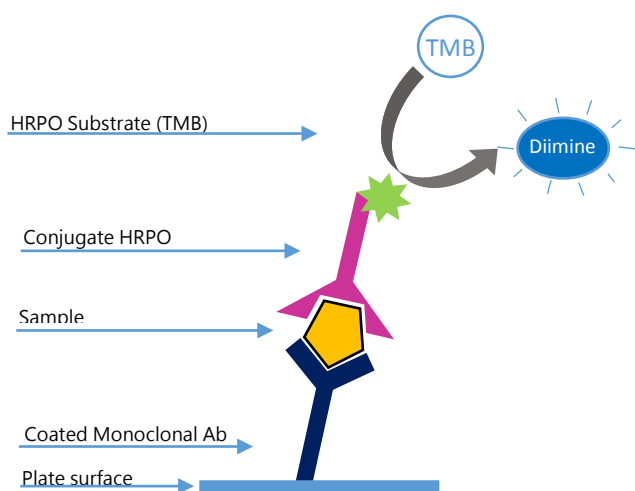INGEZIM PCV DAS is an enzymatic assay based on a Double Antibody Sandwich ELISA technique, which uses two monoclonal antibodies (MAb) specific of VP2 of PCV2. One for coating and another one as conjugate.

TECHNICAL BASIS OF THE KIT

1. Plates are coated with a MAb specific to PCV's nucleoprotein.
2. If the samples contain protein N of PCV, the MAb will capture this protein.
3. When another MAb-PO, specific to VP2 of PCV is added, it will bind to the protein. The binding is detected by the development of a colorimetric reaction after the addition of the substrate.

APPLICATION

Control of the viral load in vaccines by titration.

INTERPRETATION OF THE RESULTS

The titre of the sample is established as the last dilution with an OD higher than the negative control OD multiplied by a fixed factor.

VALIDATION

SENSITIVITY

The assay is able to detect **3ng** of purified recombinant VP2 of PCV type II.

ESPECIFICITY

The assay has been evaluated against different recombinant proteins of other viruses not related to PCV (VP7 of BTV4, BTV8, BTV1, ADV gB, CCV, p80 of BVD), foetal calf serum, Leukocytes and plasma. No cross-reaction was found.

COMPOSITION OF THE KIT

- Microtitration plates of 96 wells
- Vials with Positive Control
- Vials with Negative Control
- Vials with Peroxidase Conjugate
- Bottle with Washing Solution
- Bottle with diluent
- Bottle with stop solution
- Bottle with substrate (TMB) ready to use

PRODUCT MANUFACTURED BY INGENASA

SHELF LIFE: 12 months
Stored at 2-8°C

Ed.120117